

updater.zip

Contains files used to update the device and the PC software.

Procedure:

1) Unzip the zip file. You get a folder "updater" and a program file "trace_mon.exe". The updater is designed to update the unit and the program file is the respective software for the control PC.

2) Copy the folder "updater" in the root directory of a USB stick

3). Insert the USB stick into the TRACE C2 Control and turn on the device. After booting (white LED), the LED changes to red. Turn it off again and repeat the process until after booting a traffic light signal red, yellow, green appears.

(Note: The traffic signal serves as verification that the USB stick has been detected properly. Sometimes USB memory sticks boot too slowly and it may happen that the TRACE C2 Control does not realize it has a USB Stick with further instructions on board. In this case several attempts are necessary.)

4) To update the program file overwrite the existing file "trace_mon.exe" with the new version in your program directory on your PC.